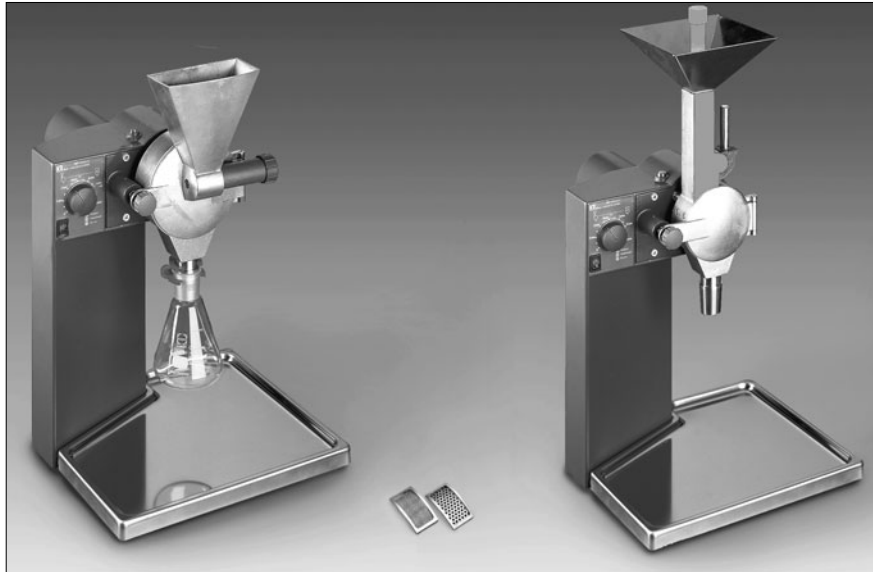
# IKA MF 10

<i>BETRIEBSANLEITUNG</i>	<i>D</i>	<i>6</i>
<i>OPERATING INSTRUCTIONS</i>	<i>GB</i>	<i>16</i>
<i>MODE D'EMPLOI</i>	<i>F</i>	<i>26</i>
<i>INSTRUCCIONES DE MANEJO</i>	<i>E</i>	<i>36</i>
<i>HANDLEIDING</i>	<i>NL</i>	<i>46</i>
<i>ISTRUZIONI PER L'USO</i>	<i>I</i>	<i>56</i>
<i>DRIFTSANVISNING</i>	<i>S</i>	<i>66</i>
<i>DRIFTSINSTRUKS</i>	<i>DK</i>	<i>76</i>
<i>DRIFTSVEJLEDNING</i>	<i>N</i>	<i>86</i>
<i>KÄYTTÖOHJE</i>	<i>SF</i>	<i>96</i>
<i>INSTRUÇÕES DE SERVIÇO</i>	<i>P</i>	<i>106</i>

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**Insert sheet for MF 10 Operating Manual**

The following changes have been made:

**The locking bolt item A15 is eliminated. Its funktion is replaced by the socket wrench, included with delivery.**

The locking of the shaft described in the chapter on Commissioning, sub-section Impact mill / assembly of the impact grinder head or Sub-section cutting mill / assembly of the cutting grinder head is replaced by the bolt, item A15. While the cylinder screw is being tightened, the rotors are kept in place by the socket wrench, included with delivery. To do this, the wrench is applied so that the sides of the wrench reach in and grip inside the front side drill holes of the rotors. For the hammer rotor, the disk must be rotated so that the drill holes in the rotor are accessible for the sides of the wrench.

**The feather key, item A7, is permanently installed in the rotor shaft.**

The assembly or disassembly of the feather groove, described several times in the chapter on commissioning, is eliminated. For installation or disassembly of the heads, the user must take care that the feather groove is in the area of the cut-out (on the rear side of the grinding track). Then it will be possible to pull the head off forwards or to set it back in place.

**Spare parts list:** MF10 basic: Items 15; 16; 17 and 49 are eliminated.

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## Contents

	Page
Safety instructions	16
Unpacking	16
Important information	17
Proper use	17
Commissioning	18
Troubleshooting	22
Maintenance and cleaning	23
Associated standards and regulations	24
Accessories	24
Garantie	24
Technical data	25
List of spare parts MF 10 basic / Gearing	116
Spare parts diagram MF 10 basic	117
Spare parts diagram Gearing	118
Spare parts diagram MF 10.1	119
List of spare parts MF 10.1 nd. MF 10.2	120
Spare parts diagram MF 10.2	121

## Safety instructions

- The machine may only be opened by trained specialists - even during repairs. The machine is to be unplugged from the mains before opening. Live parts inside the machine may still be live for some time after unplugging from the mains.  
Make certain that the mill drive stands firmly on a table.
- The mill is to be operated only under the supervision of an authorized operator.
- The mill is to be switched off completely using the rocker switch (Pos. A13) before opening the door.
- Excess pressure develops with closed collecting vessels, in particu-

lar when using dry ice or liquid nitrogen. This can lead to ground material escaping upwards out of the funnel.

- Operating personnel must wear protective personal equipment during the grinding procedure as fragments from the material to be ground may escape.
- Feeding stock residue that remain in the funnel may not be manually removed during operation or with unsuitable tools e.g. screwdrivers. They should only be removed with suitable cleaning tools when the device is switched off.
- Fingers and/or other articles should not be placed into the discharge cone (Pos. P1/S1) in order to clean it during operation.
- Cleaning is only permitted when the device is switched off and the power supply plug has been disconnected.
- No highly combustible materials may be cut up.

**Note:** After longer grinding cycles the grinding chamber and the crushing tools may be hot. Let the mill cool off before cleaning or performing any repair work.

**Note:** Exercise extreme caution when assembling, disassembling or cleaning the cutting-grinding head, particularly with regard to the sharp blade edges on the rotor (Pos. S20). These parts are to be handled with extreme care.

Wear protective personal equipment.

## Unpacking

Please unpack the device carefully and check it for damage. It is important that any transport damage is detected when the device is unpacked. If necessary, an inventory of the damage should be made immediately (postal service, railway, haulage company).

The delivery quantity for the drive includes: an MF10 drive unit, a collecting pan, a mains cable, a feather key, an Allen wrench, a fork wrench and an operator's manual.

The delivery quantity for the impact grinding head includes: an impact grinding surface with built-on door and three screwed-in fastening pins, a dosing sluice, a hammer rotor, a countersunk screw, a disk, a cover and

the filter which you ordered.

The delivery quantity for the cutting-grinding head includes: a cutting-grinding head with three built-in bed knives, built-on door, tabular slide valve and funnel, as well as three screwed-in attaching pins, a plunger, a blade rotor with three mounted rotor blades, a cheese head screw, a disk and the filter which you ordered.

## Important information

The MF 10 analytical mill is a continuously operating universal mill. It makes it possible to use two different grinding methods which cover the greatest range of all grinding tasks occurring in the laboratory.

### **Impact grinding:**

Application for hard, brittle and/or dried materials or ground materials embrittled as a result of cold (e.g. grain, coffee, coal, etc.).

In this case, the ground material is cut up by means of rotary beaters and comminuted on the grinding surface attached to the blow bars. The ground material is broken and ejected into a suitable collecting pan by means of corresponding filters. The feed dosage provides for optimal dosage so that the mill can work at the optimal operating point. The grinding chamber is locked so that it is dustproof, the collecting pans are attached with standard ground joints. Thus the loss ground material is kept to a minimum.

### **Cutting-Grinding:**

Application for soft fibrous ground material (e.g. hay, straw, plastic waste, paper, wood, etc.).

In this case, the ground material is cut between rotary blades and standing blades attached at the grinding surface. The ground material is conveyed into the collecting pan by means of filters here as well. The feed unit (guide channel) makes it possible to load awkward, bulky material. The cutting loads mean that less energy is introduced into the ground material and thus results in less heat generation. This keeps ground material loss to a minimum.

## Proper use

The area of application for the universal mill is very extensive, the list of materials which are capable of being ground and indicated here is not exhaustive. The scope of application can be extended further by grinding attempts with the respective grinding method and setting the appropriate grinding parameters (rotor speed, filter hole diameter, sample preparation e.g. cooling).

### **Impact grinding:**

The hammer mill crushes soft, medium-hard and brittle materials to approximately. Mohs hardness 6 (feldspar: 6 Mohs, quartz: 7 Mohs). With greater hardness the wear on the grinding tools is much higher and grinding is therefore uneconomical. Anything can be ground that breaks, is dry does not contain very fatty materials.

The following is a selection of substances drying to grind: grain, barley, corn, malt, pectin, roasted coffee, nutshells, bones, ergot, coal, charcoal, coke, peat, ash, cellulose, drugs, artificial fertilizers, refractory clay, feed, spices, glass, resin, potash, cores, salts, cinders, stones and soil (bauxite, gypsum, limestone, dry clay, mica, graphite, pumice stone, gneiss, magnetite, heavy spar, sandstone, brown iron ore, etc.)

The scope of application can be further enhanced through embrittlement of the ground material (e.g. with dry ice and/or liquid nitrogen). Tough ground material must be cooled, e.g. crushed dry ice put into the grinding funnel.

The ground material must be coarse-crushed into maximum 10mm chunks beforehand.

**Note:** With abrasive materials (e.g. stones or glass) it will be necessary to clean the dosing sluice repeatedly because settles grinding dust in the guides and it thus rotating the dosing sluice is no longer is possible.

### **Cutting-Grinding:**

The cutting-grinding mill cuts up bulky, elastic, fibrous, cellulose and soft materials. Mixing materials such as various types of waste must be free of iron and non-ferrous metals. The feedings may not be damp and greasy, as they will otherwise stick to the grinding surface and filter.

The following is a list of the materials that may be processed: Leaves, fibers, spices, grain, hop, wood, pasteboard, paper, hay, straw, plastics, tobacco, peat, feed, pasta, roots and branches, etc.

The scope of application can be further enhanced through embrittlement of the ground material (e.g. with dry ice and/or liquid nitrogen).

Bulky material is metered with the tubular slide valve, the feed grain size should not be not greater than 15mm. Bulky materials (e.g. plastic bags) are metered by means of plungers, but the pieces should not be too large (e.g. pieces of fir wood: max. diameter 10mm, max. length 40mm; pieces of foil: max. width 30mm, max. length 100mm).

## Commissioning

Make certain that the mill drive stands firmly on a table. Comply with the ambient conditions given in the technical data.

The collecting pan (Pos. A19) is placed on its base in such a way that the round magnets fit into the recess on the base.

### **Impact grinding mill:**

#### ***Assembling the impact grinding head:***

The ON/OFF switch Pos. A13 must be set to zero and the power supply plugs must have been disconnected out before beginning assembly of the impact grinding head.

First the three threaded bolts Pos. P5 are screwed out to the same height (distance between hexagon of the bolt and the flat surface on the grinding head is approximately 2mm, this corresponds to two turns).

Make sure that the flange piece on the drive and on the head are clean and free of dust. Abrading both fitting parts with a slightly oiled rag facilitates assembly. The head is now pushed onto the drive, while making certain that the parts do not jam. The threaded bolt ends must fit into the drillings of the transmission cover. Thus the head has been rotated slightly in a clockwise direction. After the bolts are fitted in the head is turned until it is stopped by the bolts and so that the head now stands upright.

Now close and lock the grinding chamber door with the operating knob Pos. P10 (3 to 5 revolutions are sufficient). This ensures that the grinding chamber is optimally aligned so that the door can be easily opened and closed later.

Now the screw mountings are tightened with the fork wrench provided. Make certain that the pins are tightened toward the head.

**Note:** **Tighten the pins evenly** (see Fig. 1)!

Now the hammer rotor Pos. P13 can be installed. For this the fitting key Pos. A7 is first inserted into the groove of the drive shaft Pos. G6, make certain that the groove is free from dirt and dust here as well. The rotor is pushed onto the drive shaft as shown in Fig. 3, fastened with the disk Pos. P11 and the countersunk screw Pos. P15. The countersunk screw is tightened with the Allen wrench provided, while the pin Pos. A15 is pressed. The pressed pin prevents the shaft from slipping while the countersunk screw is tightened. The pin engages in two positions. These can be found by pressing the pin and at the same time turning the shaft (simply turn the rotor with the tool until the pin engages).

**Note:**

After tightening the cheese head screw pin Pos. A15 release and turn the hammer rotor once (360°) using the Allen wrench. The rotor may not touch. Repeat assembly if necessary.

Now the filter Pos. Z1 that is suitable for the grinding procedure can be inserted; to do so the filter is pushed into the grinding chamber from the front. The filter and the filter holder surface in the grinding chamber must be free from dirt and dust.

Now the grinding chamber door can be closed and locked with the operating knob Pos. P10. The screw mounting of the operating knob is screwed in by pressing and twisting it into the internal thread on the drive cover. The first turns are easy, if turning becomes difficult then the door may be closed, but the safety switch is not yet unlocked. Continue turning until stop (approximately two revolutions), only then is the safety switch unlocked and the mill ready to start.

### **Grinding operation with impact grinding mill:**

The suitable filter is pushed in and the door is locked before beginning the grinding procedure.

Afterwards a collecting pan is fastened to the discharge cone (standard ground joint NS 29). Suitable containers are e.g. Erlenmeyer flasks Pos. Z3, round-bottomed flasks, test tubes made of either glass or plastic. The containers are to be secured with a fixing clip Pos. Z2.

For heat sensitive materials or with ground materials that have a very high increase in volume during cut up, it is advisable to use a two-necked flask with a filter attached to its second (e.g. cotton wool, etc...). The ground material is cooled from the air flow and no excess pressure is created. For very critical ground materials (very fine powder after being cut up) the ground material can also be sucked off by means of a commercial vacuum cleaner with a suction tube adapted at the discharge cone. The powder is then in the vacuum cleaner bag where it can be removed for further processing steps.

The ground material which has prepared to the correct feed grain size is filled into the funnel and then the cover Pos. P17 is applied. The speed knob Pos. A12 is turned to the left stop (3000 1/min).

The mill is started by shifting the rocker switch Pos. A13 to Position 1. First the green „Power“ LED is lit and then „Overload“ LED lights up briefly before the motor smoothly starts. The speed can be set by turning the speed knob to the desired speed. A speed of more than 4500 1/min is recommended for impact grinding.

Now the ground material can be metered by turning the dosing sluice Pos. P6 in a clockwise direction. You can hear if ground materials gets into the grinding chamber and if too much materials has been metered then the rotor speed decreases. This drop in speed is clearly indicated by the lower grinding noises and the orange „Overload“ LED lights up. Metering of the ground material should only be resumed when the speed has returned to the adjusted value or if the „Overload“-LED is no longer lit.

If this is done then the mill becomes overloaded and switches off after some time because of overheating. This is then indicated by the red „Error“ LED. In addition, a beep sounds to indicate that the mill should be switched off using the rocker switch Pos. A13. The mill must then

cool off for a period in order to be able to start in again (see chapter on Troubleshooting).

In addition, the ground material is no longer cut up the lower speed when the unit is operated with an overload because the impact energy no longer suffices for breaking the ground material. Keeping the ground material in the grinding chamber and the ground material to heat up. In extreme cases this can result in the ground material caking on the grinding chamber parts which then entails a time-consuming cleaning operation.

If the funnel is empty then the mill remains turned until all of the ground material from the grinding chamber is in the container. This is the case if cutting up noises can no longer be heard.

When it comes to cleaning the grinding chamber experience has shown that it makes sense to briefly switch the unit to maximum speed after the grinding operation has been completed and only then to turn the speed knob to the left stop. The mill is then shut off with the switch.

The filter is removed after opening the grinding chamber door and any remaining ground material is wiped into the collecting pan.

**Note:** After longer grinding cycles the grinding chamber and the filters may have an increased temperature. Protective personal equipment should be worn in this case.

It may help to crush the ground material in several steps, i.e. a filter with large perforations is used for pre-crushing and finer ones in each of the following steps.

Advantages:

- fewer failures, e.g. due to clogged filters or overloading the motor
- less heat-up of ground material since the retention time of the ground material in the grinding chamber is substantially reduced
- closer grain size spectrum for the final material
- normally less total grinding with better results as despite several runs comminution is performed more quickly and more effectively

### ***Disassembling the impact grinding head:***

Before disassembling the head of the mill, the mill itself must first cool down as the grinding chamber and the beater rotor cannot be detached otherwise.

The power supply plug must be disconnected before disassembly. After the door is opened the countersunk screw Pos. P15 is loosened by means of an Allen wrench. For this the shaft is blocked by pressing the pin Pos. A15. If the hexagon socket is dirty then it must be cleaned by means of a brush. After the countersunk screw and the disk Pos. P11 are removed the hammer rotor Pos. P13 is taken off and the fitting key Pos. A7 removed. Now the three screw mountings Pos. P5 are loosened by means of a fork wrench and the grinding chamber is rotated in a clockwise direction and taken off.

### **Cutting mill:**

#### ***Assembling the cutting mill head:***

**Caution:** The blades on the rotor and in the grinding chamber have sharp edges. Both parts should only be handled when wearing protective gloves.

The ON/OFF switch Pos. A13 must be zero and the power supply plugs must have been disconnected out before beginning assembly of the cutting-grinding head.

First the three threaded bolts Pos. S12 are screwed out to the same height (distance between hexagon of the bolt and the flat surface on the grinding head is approximately 2mm, this corresponds to two turns).

Make sure that the flange piece on the drive and on the head are clean and free of dust. Abrading both fitting parts with a slightly oiled rag facilitates assembly. The head is now pushed onto the drive, while making certain that the parts do not jam. The threaded bolt ends must fit into the drillings of the transmission cover. Thus the head has been rotated slightly in a clockwise direction. After the bolts are fitted in the head is turned until it is stopped by the bolts and so that the head now stands upright.

Now close and lock the grinding chamber door with the operating knob Pos. S9 (3 to 5 revolutions are sufficient). This ensures that the

grinding chamber is optimally aligned so that the door can be easily opened and closed later.

Now the screw mountings are tightened with the fork wrench provided. Make certain that the pins are tightened toward the head.

**NOTE:** **Tighten the pins evenly** (see Fig. 1) !

Now the hammer rotor Pos. S20 can be installed. For this the fitting key Pos. A7 is first inserted into the groove of the drive shaft Pos. G6, make certain that the groove is free from dirt and dust here as well. The rotor is pushed onto the drive shaft as shown in Fig. 2, fastened with the disk Pos. S22 and the cheese head screw Pos. S21. The countersunk screw is tightened with the Allen wrench provided, while the pin Pos. A15 is pressed. The pressed pin prevents the shaft from slipping while the countersunk screw is tightened. The pin engages in two positions. These can be found by pressing the pin and at the same time turning the shaft (simply turn the rotor with the tool until the pin engages).

**Note:**

After tightening the cheese head screw pin Pos. A15 release and turn the hammer rotor once (360°) using the Allen wrench. The rotor may not touch. Repeat assembly if necessary.

Now the filter Pos. Z1 that is suitable for the grinding procedure can be inserted; to do so the filter is pushed into the grinding chamber from the front. The filter and the filter holder surface in the grinding chamber must be free from dirt and dust.

Now the grinding chamber door can be closed and locked with the operating knob Pos. S9. The screw mounting of the operating knob is screwed in by pressing and twisting it into the internal thread on the drive cover. The first turns are easy, if turning becomes difficult then the door may be closed, but the safety switch is not yet unlocked. Continue turning until stop (approximately two revolutions), only then is the safety switch unlocked and the mill ready to start.

### **Grinding operation with the cutting mill:**

The suitable filter is pushed in and the door is locked before beginning the grinding procedure.

Afterwards a collecting pan is fastened to the discharge cone Pos. S1 (standard ground joint NS 29). Suitable containers are e.g. Erlenmeyer flasks Pos. Z3, round-bottomed flasks, test tubes made of either glass or plastic.

The containers are to be secured with a fixing clip Pos. Z2.

For heat sensitive materials or with ground materials that have a very high increase in volume during cut up, it is advisable to use a two-necked flask with filter attached to its second (e.g. cotton wool, etc.). The ground material is cooled from the air flow and no excess pressure is created. For very critical ground materials (very fine powder after being cut up) the ground material can also be sucked off by means of a commercial vacuum cleaner with a suction tube adapted at the discharge cone. The powder is then in the vacuum cleaner bag where it can be removed for further processing steps.

The speed knob Pos. A12 is turned to the left stop (3000 1/min).

The mill is started by shifting the rocker switch Pos. A13 to Position 1. First the green „Power“ LED is lit and then „Overload“ LED lights up briefly before the motor smoothly starts.

The speed can be set by turning the speed knob to the desired speed. A speed up to 5000 1/min is recommended for cutting-grinding while higher speeds should only be selected in exceptional cases (e.g. greater energy requirements such as for dry pieces of wood). The optimal speed must be determined by way of experiment. The higher the speed, the greater the cutting-grinding power. However, more energy is introduced into the ground material and thus the thermal load is increased.

The plunger Pos. S6 has to be removed from the guide channel before the ground material is fed in through the funnel Pos. S5. Make certain that the tubular side valve Pos. S8 is closed. Bulky materials can be simply filled into the funnel.

Metering is carried out by lifting the tubular slide valve.

Bulky material should be fed into the guide channel with the tubular side valve closed and the plunger must be fitted to in front of the tubular side valve in order to prevent chunks of ground material from spurting out.

The plunger may only be pulled out when the motor is running if the tubular side valve is closed.

**WARNING: Wear protective personal equipment (protective glass, etc.)**

You can hear if ground material gets into the grinding chamber and if too much material has been metered then the rotor speed decreases. This drop in speed is clearly indicated by the lower grinding noises and the orange „Overload“ LED lights up. Metering of the ground material should only be resumed when the speed has returned to the adjusted value or if the „Overload“ LED is no longer lit.

If this done then the mill becomes overloaded and switches off after some time because of overheating. This is then indicated by the red „Error“ LED. In addition, a beep sounds to indicate That the mill should be switched off using the rocker switch Pos. A13. The mill must then cool off for a period in order to be able to start it again (see chapter on Troubleshooting).

When it comes to cleaning the grinding chamber experience has shown that it makes sense to insert the plunger into the guide channel as far as possible and then to briefly switch the unit to maximum speed after the actual grinding operation has been completed and only then to turn the speed knob back to the left stop. The mill is then shut off with the switch.

Bulky and fibrous ground material, e.g. straw, hay, etc. should not be fed into the funnel lengthwise or in the form of larger positions. The plunger should be used to introduce them into the grinding chamber. Note that the plunger is coded and thus only fits into the guide channel in one direction (edge bevelling) on the plunger and in the guide channel have to match each other). The „V“ on the plunger must be visible from the front. It has proved to be advantageous in some cases

if the plunger is turned around and the bulky material is pushed from the funnel into the guide channel with the cylindrical end of the plunger.

The better a sample is prepared (pre-crushed to the optimal feed size), the easier the metering and the faster the grinding operation. In addition, failures such as rotor blocking can be virtually eliminated.

The filter is removed after opening the grinding chamber door and any remaining ground material is wiped into the collecting pan.

**Note:**

After longer grinding cycles the grinding chamber and the filter s may have an increased temperature. In this casewait until the parts have cooled off.

It may help to crash the ground material in several steps, i.e. with large perforations is used for pre-crushing and finer ones in each of the following steps.

Advantages:

- fewer failures, e.g. due to clogged filters or overloading the motor
- less heat-up of ground material since the retention time of the ground material in the grinding chamberchamber is substabtily reduced
- closer grain size spectrum for the final material
- normally less total grinding with better results as despite several runs comminution is performed more quickly and more effectively

**Disassembling the cutting-grinding head:**

**Warning:**

The blades on the rotor and in the grinding chamber have sharp edges. Observe safety regulations.

Before disassembling the head of the mill, the mill itself must first cool down as the grinding chamber and the beater rotor cannot be detached otherwise.

The power supply plug must be disconnected before disassembly.

After door is openend the countersunk screw Pos. S21 is loosened by means of Allen wrench. For this the shaft is blocked by pressing the pin Pos. A15. If the hexagon socket of the screw is dirty then it must be cleaned by means of a brush. After the cheese head screw and the disk Pos. S22 are removed the hammer rotor Pos. S4 is taken off and the fitting key Pos. A7 removed. Now the three screw mountings Pos. 12 are loosened by means of a fork wrench and the grinding chamber is rotated in a clockwise direction and taken off.

## Troubleshooting

**Mill does not start:**

Possible causes:

Power supply plug is not correctly plugged into the equipment and/or the socket („Power“ LED does not light up if the rocker switch Pos. A13 is moved to Position 1).

Grinding Chamber door not properly locked, („Power“ LED, „Error“ LED are lit and beep is heard when the rocker switch Pos. A13 is moved to Position 1).

Remedy:

First shift the rocker switch Pos. A13 to 0. Check voltage supply and the power supply plug is necessary.

Or if „Error“ LED is lit and beeping can be heard close the grinding chamber door and tighten the turning handle Pos. P10 / S9 until stop.

The rocker switch should only then be shifted to Position 1 again.

**Engine hums after the rocket switch Pos. A13 is shifted but the rotor does not begin to run.**

Possible causes:

Ground material chunks is in the grinding chamber before grinding begins and blocks the rotor.

Switch the mill off at the rockers switch Pos. A13 and disconnect the power supply plug. Open the door and remove the ground material chunks from the grinding chamber. After the door is closed and the mill is connected to the mains again, you can switch the mill on again using the rocker switch Pos. A13.

**Engine switches off during operation:**

Possible causes:

Overloading of the engine because of too large feeding amounts which leads to heating up of the engine. Thus the motor thermal protective switch is triggered and switches the mill off. The overheating of the motor is indicated by the red „Error“ LED becoming lit and a beep is heard as well.

Rotor blocking as a result of feed chunks that are too large or too tough.

Rotor blocking due to improper speed adjustment (in general, the speed is too low for the given application).

Remedy:

Switch the mill off at the rocker switch Pos. A13 and disconnect the power supply plug. Open the door and remove any blocked ground material or ground material that has been cut up.

If the disturbance occurred because of overloading, let the mill cool off before you start it again.

## Maintenance and cleaning

The grinding chamber interior (funnels, dosing equipment, rotor, grinding surface, filter and outlet nozzle) with a brush (no wire brush) should be cleaned after the grinding procedure. You can also use rags or cleaning brushes. The grinding surface does not need to be disassembled for cleaning between grinding operations with the same feed material.

If a new grinding procedure with other feedings is to be performed whereby no contamination may be produced by the previously ground material, then more thorough cleaning is necessary which means dismantling the grinding chamber and the rotors. See the chapter „Commissioning“ for assembly and disassembly.

The grinding chamber parts can be washed with water and, if required, with a conventional dish-washing agent.

Be certain that the parts are dry before being reassembled.

**Caution:**

Refer to the chapter on „Safety Instructions“ when cleaning the cutting-grinding surface and the blade rotor as the screwed on blades are very sharp and may result in injuries if they are improperly handled.

Clogged filters can be cleaned with brushes or compressed air. If ground material parts have settled into the filter holes in such a way that cleaning alternatives are not sufficient, then you will have to clean the clogged holes with suitable needles or clean the filter in an ultrasonic bath.

If beaters or blades and the grinding surface become worn in the course of the time, particularly with constant grinding of very hard substances, so that quick operation with the appropriate quality is no longer possible, then the worn out parts should be replaced.

**Blades may not be reground as the blade clearance will become too large.**

**Fore reasons of safety, only new, original IKA bolts may be used for fastening the knives.**

The mill IKA MF 10 is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate.

**When ordering spare parts, please give the manufacturing number shown on the type plate, the machine type and the name of the spare part.**

Please send in equipment for repair only after it has been cleaned and is free from any materials which may constitute a health hazard. Use only cleansing agents which have been approved by IKA to clean IKA devices. To remove use:

Dyes	isopropyl alcohol
Construction materials	water containing tenside / isopropyl alcohol
Cosmetics	water containing tenside / isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

For materials which are not listed, please request information from IKA. Wear the proper protective gloves during cleaning of the devices.

Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does not destroy the instrument.

## Associated standards and regulations

### Associated EU guidelines

EMC-guideline: 89/336/EWG  
Low voltage-guideline: 73/023/EWG  
Machine-guideline: 89/392/EWG

### Construction in accordance with the following safety standards

EN 61 010-1	EN 61 010-2-51
EN 292-1,	EN 292-2
EN 60 204-1	EN 414
UL 3101-1	CAN/CSA C22,2 (1010-1)

### Construction in accordance with the following EMV standards

EN 55 014-1	
EN 50 081	EN 50 082-1

## Accessories

### Sieve insert with hole sizes:

Ø 0,25mm	Ø 1,50mm
Ø 0,50mm	Ø 2,00mm
Ø 1,00mm	Ø 3,00mm

<b>MF 10.1</b>	Cutting mill head
<b>MF 10.2</b>	Impact grinding head

## Guarantee

You have purchased an original IKA laboratory machine which meets the highest engineering and quality standards.

In accordance with IKA guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

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### MF 10 basic

Item Designation

2	Graphit packing	20	Device base
3	BLP logic	22	Base
4	Device plug	28	Ring
5	Housing cover	29	Pulse recorder
6	BLP output	30	Round wire snap ring
7	Fitting key	31	O-Ring
8	Cap	32	Indented ring
11	Front foil	35	Triac
12	Operating knob	36	Square disk
13	Rocker switch	37	Ribbon cable
15	Bolt	49	O-Ring
16	Pressure spring	50	Clip
17	O-ring	51	Temperature switch
18	Housing front part	2001	Gearing; complete
19	Collecting pan	3001	Fork-type light barrier

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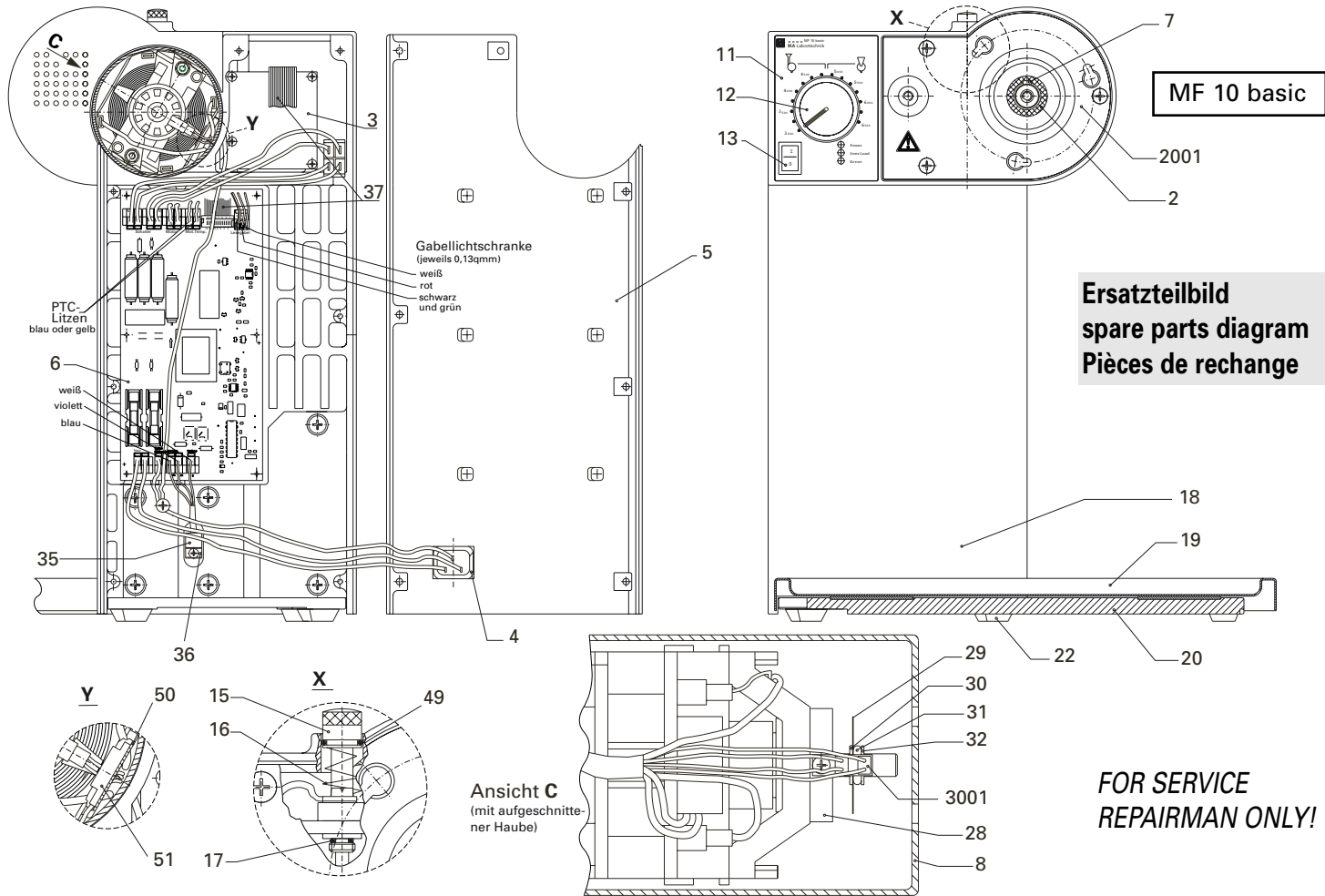
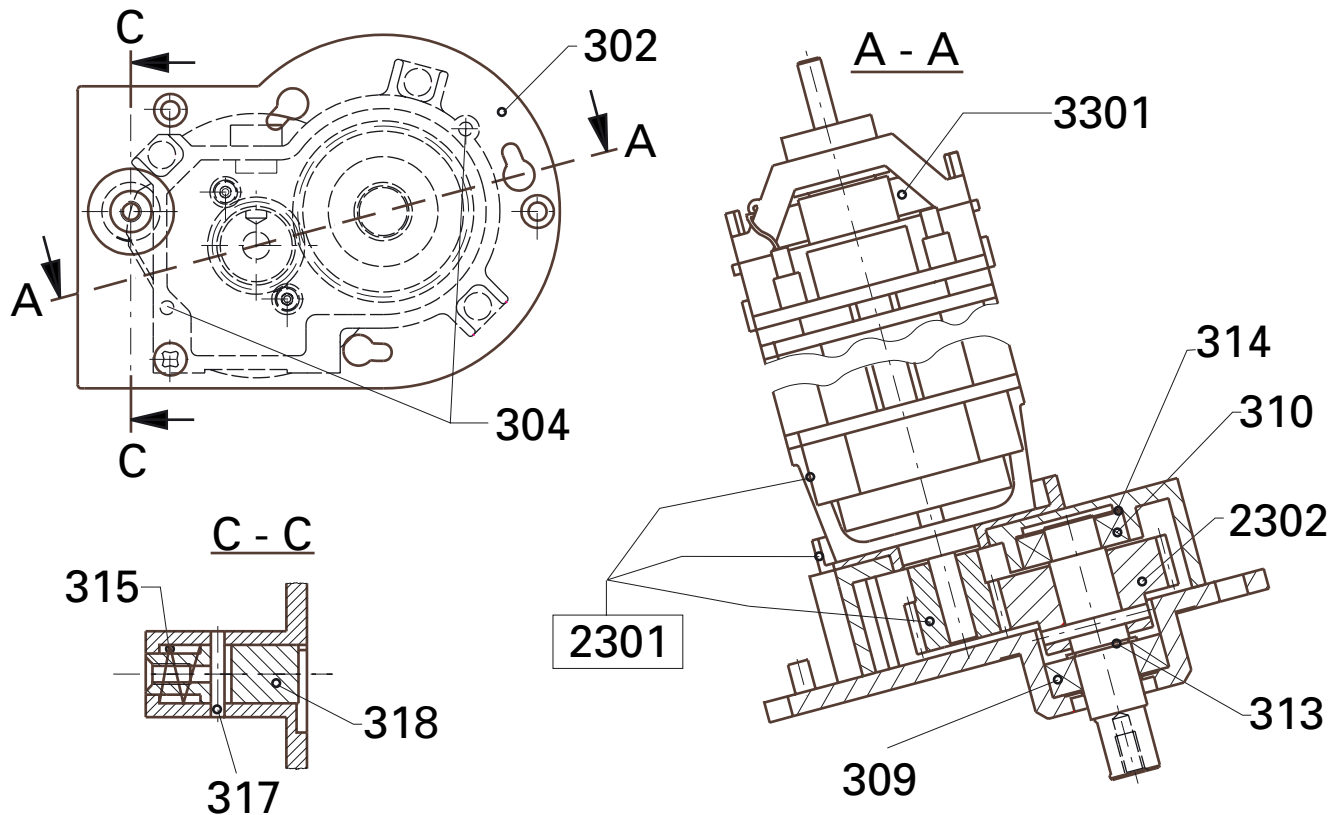
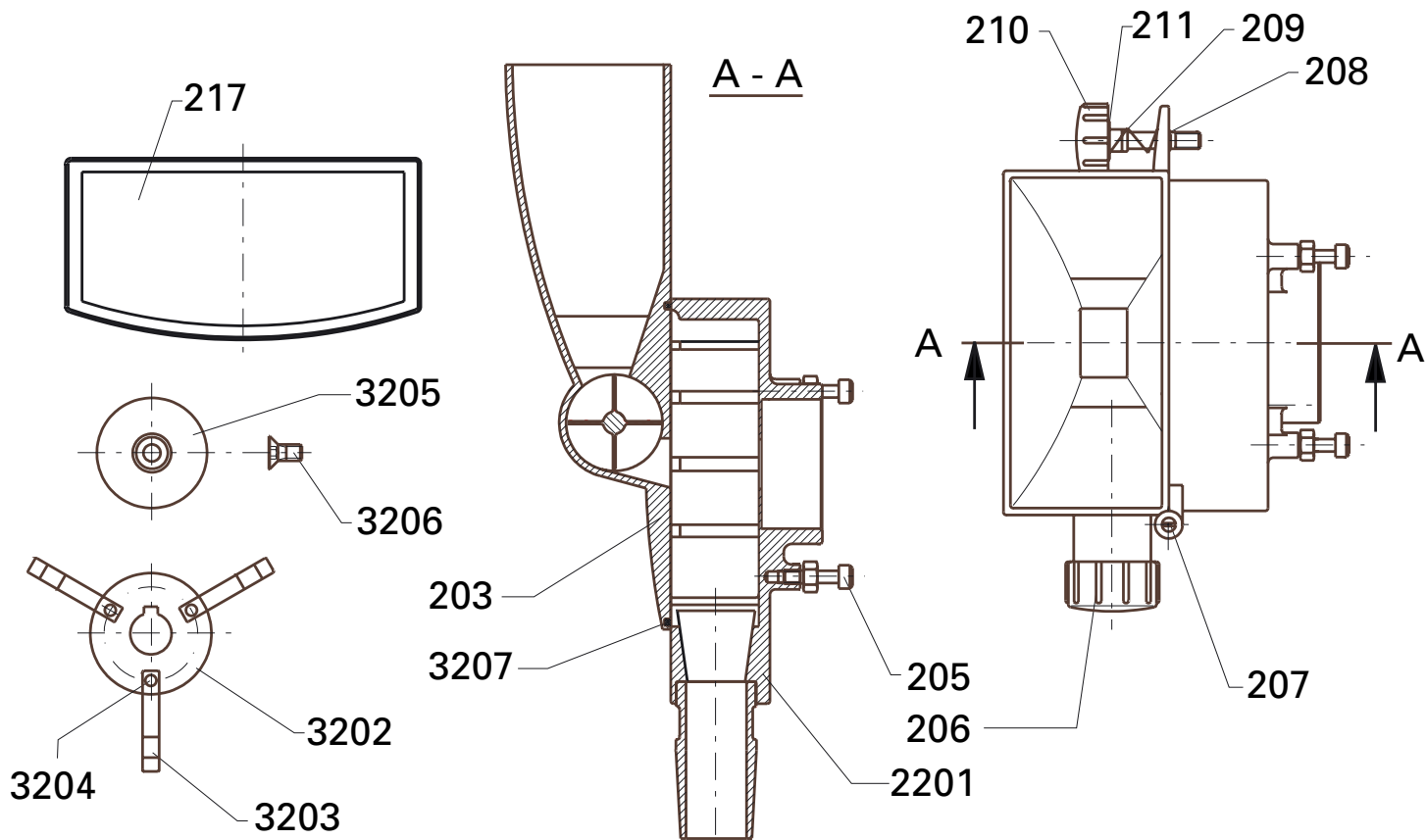
## **IKA® Works, Inc.**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

### North America

---

**IKA® Works, Inc.**  
2635 North Chase Pkwy SE  
Wilmington, NC 28405-7419  
USA  
Tel.: 800 733-3037  
Tel.: +1 910 452-7059  
Fax: +1 910 452-7693  
E-mail: usa@ika.net  
www.ika.net

## **IKA® Works (Asia) Sdn Bhd**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

### Asia - Australia

---

**IKA® Works (Asia) Sdn Bhd**  
No. 17 & 19, Jalan PJU 3/50  
Sunway Damansara Technology Park  
47810 Petaling Jaya  
Selangor, Malaysia  
Tel.: +60 3 7804-3322  
Fax: +60 3 7804-8940  
E-mail: sales@ika.com.my  
www.ika.net

## **IKA® Japan K.K.**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

### Japan

---

**IKA® Japan K.K.**  
293-1 Kobayashi-cho  
Yamato Koriyama Shi, Nara  
639-1026 Japan  
Tel.: +81 743 58-4611  
Fax: +81 743 58-4612  
E-mail: info@ika.ne.jp  
www.ika.net

## **IKA® Works Guangzhou**

LABORATORY TECHNOLOGY  
ANALYZING TECHNOLOGY  
PROCESSING EQUIPMENT

### China

---

**IKA® Works Guangzhou**  
173 - 175 Friendship Road  
Guangzhou Economic and Technological  
Development Zone  
510730 Guangzhou, China  
Tel.: +86 20 8222-6771  
Fax: +86 20 8222-6776  
E-mail: sales@ikagz.com.cn  
www.ika.net