MR Hei-Mix L

Betriebsanleitung
Instruction Manual
Mode d'Emploi
Instrucciones de
Empleo
Istruzioni per l'uso

D	DEUTSCH	Seite	3 - 12
E	ENGLISH	page	13 - 22
F	FRANCAISE	page	23 - 32
ES	ESPAÑOL	página	33 - 42
I	ITALIANO	pagina	43 - 52

INHALT

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+ EN 61010-2-051

SUMMARY

SCOPE OF DELIVERY	14
GENERAL INFORMATION	14
SAFETY INFORMATION	15
INTENDED USE	16
SET-UP	16
1 Unpacking	16
2 Setting up the instrument	16
3 Electrical connection	16
OPERATION	17
4 Switching the instrument on	17
5 Setting the stirring speed	17
CLEANING AND MAINTENANCE	17
TRANSPORT AND STORAGE	18
DISPOSAL	18
TROUBLESHOOTING	18
REPLACEMENT PARTS AND ACCESSORIES.....	19
SPECIFICATIONS.....	19
WARRANTY, LIABILITY & COPYRIGHT	20
QUESTIONS, REPAIRS.....	21
APPLICABLE STANDARDS AND DIRECTIVES	22

SCOPE OF DELIVERY

Product	Quantity	P/N	
		230 V / 50/60 Hz	115 V / 60 Hz
<i>Hei-Mix L</i>	1	505-00000-00	505-00000-01
Instruction manual	1	01-005-004-65	01-005-004-65
AC power cord	1	14-007-003-81	14-007-003-89

GENERAL INFORMATION

Read the instruction manual with care and ensure that all users read it with care prior to using the instrument.

Please store the instruction manual in a place easily accessible to every user.

A EURO AC plug (DIN 49441 CEE 7/VII 10/ 16 A 250 V) is standard on all instruments.

For North America, instruments feature standard US plugs (NEMA Pub. No. WDI.1961 ASA C 73.1 . 1961 page 8 15A 125V).

When operating the instrument in countries with different AC plug systems, use an approved adapter or have a qualified electrician replace the AC plug with an approved model suitable for the country of operation.

The instrument is earthed as supplied. When replacing the original AC plug, ensure that the earth conductor is connected to the new plug.

SAFETY INFORMATION

Please comply with all safety and accident-prevention regulations applicable to laboratory work.

Use great caution when working with flammable substances. Observe the relevant safety data sheets.

When connecting the instrument to an AC power outlet, ensure that your local supply voltage matches that indicated on the instrument's rating plate.

Turn the rocker switch off when not using the instrument and before disconnecting it from its AC power outlet.

Repairs may only be performed by technicians authorised by Heidolph Instruments.

Use great caution when working in the vicinity of highly flammable or explosive substances.
The instrument is not explosion-proof.

Always connect the instrument to an earthed AC power outlet.

Caution!
Beware of the effect of the magnetic field on cardiac pacemakers and data media.

Ensure that the instrument is standing on a solid surface.

INTENDED USE

The *Hei-Mix L* is a magnetic stirrer without heating. It is designed for use in chemical and biological laboratories of industrial enterprises, universities, schools and pharmacies.

The instrument is designed to mix substances through the action of a magnetically linked stirring rod in a vessel.

To ensure maximum service life, observe the specified ambient conditions (temperature and humidity) and ensure that the instrument is not exposed to a corrosive atmosphere.

SET-UP

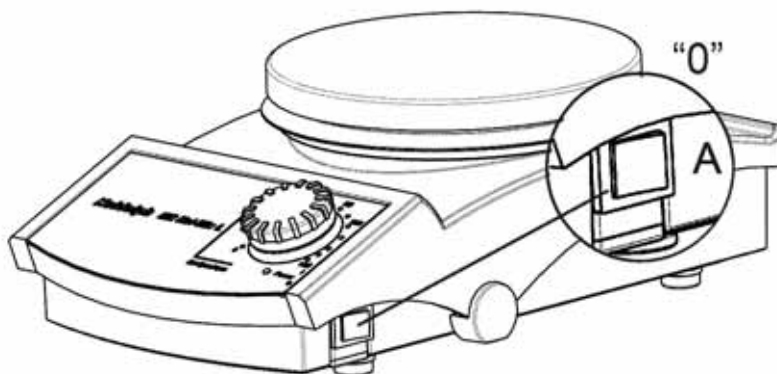
1 Unpacking

Unpack the instrument with care. Inspect for damage and report such damage or missing parts to your supplier right away.

2 Setting up the instrument

Place the instrument on a solid, horizontal surface. Keep the area around the instrument clear of other items.

3 Electrical connection

Before connecting the instrument to an AC power outlet, ensure that the rocker switch (A) is set to "0".

Plug the AC power cord into its socket (H) on the rear panel. Connect the power cord to an AC power outlet.

H

OPERATION

4 Switching the instrument on

Caution:

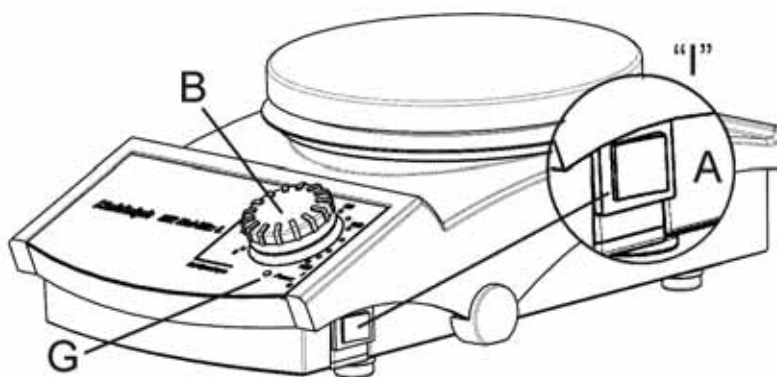
Stirring will begin immediately after the instrument is switched on in accordance with the setting of control knob (B).

Agitator will begin turning immediately!

Danger of splashing.

Before switching the unit on, turn control knob (B) fully to the left.

Turn on the instrument's power using the rocker switch (A) on the right side of the housing. The green Power LED (G) on the front panel will light.

5 Setting the stirring speed

Set the desired speed with the control knob (B) according to the scale.

The motor will start when the speed is set to 100 rpm or higher.

CLEANING AND MAINTENANCE

Use a cloth dampened with mild soapy water to clean the housing and the surface of the instrument.

Note

Do not use chlorine bleach or other chlorine-based cleaning products, abrasive cleanser, ammonia, steel wool or cleaning products with metallic components under any circumstances. These will damage the finish of the instrument.

The instrument is maintenance-free. Should repairs become necessary, please contact an authorised Heidolph Instruments repair technician. Please contact your Heidolph

Instruments dealer or a Heidolph Instruments representation for more information (see page 21).

**Caution**

Unplug the AC power cord before opening the housing.

TRANSPORT AND STORAGE

Prior to transport

Switch the instrument off and unplug the AC power cord.

Transport and storage

1. Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.
2. Store the instrument in a dry environment.

**Caution**

Do not subject the instrument to mechanical shocks or vibration while transporting it.

DISPOSAL

Please dispose of used instruments and defective components at your local recycling collection point.

Prior to disposal, sort according to materials: metal, glass, plastic, etc.

Also be sure to dispose of the packaging material in an environmentally-friendly manner.

TROUBLESHOOTING

1. Power LED (G) not lit: check power supply and AC power cord
2. No stirring function:
 - Stirring vessel does not contain a magnetic stirring rod
 - Speed control knob (B) set to "0"

REPLACEMENT PARTS AND ACCESSORIES

Accessories (optional)

Product	P/N
Heating bath, 1l - for oil up to 250 °C	504-93000-00
Heating bath, 1l, PTFE coated - for water up to 100 °C	504-93100-00
Heating bath, 2l - for oil up to 250 °C	504-92000-00
Heating bath, 2l, PTFE coated - for water up to 100 °C	504-92100-00
Heating bath, 4l - for oil up to 250 °C	504-91000-00
Heating bath, 4l, PTFE coated - for water up to 100 °C	504-91100-00
Concave block adapter for 1l round-bottom flasks	504-94000-00
Holder rod, stainless steel 1.4305 (V2A)	509-81000-00
Holder rod, stainless steel 1.4305 (V2A) (for 4l heating bath)	509-97000-00
Temperature sensor holder	509-67000-00
Magnetic stirring rod set (3 rods), cylindrical with PTFE coating	509-56000-00
MR-Hei silicone cover	23-07-06-05-59
Heating bath liquid, RT - 240 °C	515-31000-00
DrySyn MULTI complete set, 3 x 25/50/100ml inserts	505-91900-00
DrySyn MULTI-M complete set, 3 x 100/250ml inserts	505-92600-00
DrySyn triple clamp for MULTI	505-91090-00
DrySyn triple clamp for MULTI-M	505-92090-00
MR Hei holder	509-96000-00

SPECIFICATIONS

	MR Hei-Mix L
Plate material	V2A (AISI304)
Plate dimensions	145mm dia.
Speed in rpm	100 – 1.400
Speed accuracy	+/- 2%
Max. stirring quantity (water)	20l
Power consumption	20W
Permissible ambient temperature	0 - 40 °C
Permissible relative humidity	80%
Weight	2.6 kg
Dimensions (DxWxH)	277x173x94
Protection class	IP32

WARRANTY, LIABILITY & COPYRIGHT

Warranty

Heidolph Instruments provides a three-year warranty on the products described here (with the exception of consumable parts), starting from the date of shipping from the manufacturer's warehouse. This warranty covers defects in materials and workmanship.

Transit damage is excluded from this warranty.

To obtain such warranty service, contact Heidolph Instruments (phone: +49 – 9122 - 9920-68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation are not covered by this warranty promise.

Alterations to the present warranty promise need Heidolph Instruments' consent in writing.

Exclusion of liability

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

Copyright

Heidolph Instruments is the copyright holder for all texts and images in this manual.

QUESTIONS, REPAIRS

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please contact us at the following address:

For equipment repairs jobs please call Heidolph Instruments (phone: +49 – 9122 - 9920-68) or your local authorised Heidolph Instruments dealer.

**Note:**

Please return instruments to the following address only.
Returning instruments is subject to prior approval.

Heidolph Instruments GmbH & Co. KG

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Walpersdorfer Str. 12

D-91126 Schwabach, Germany

Tel.: +49 – 9122 - 9920-68

Fax: +49 – 9122 - 9920-65

E-mail: sales@heidolph.de

**Safety information**

When returning instruments for repair that have come in contact with hazardous substances, please:

- provide precise information on the relevant medium
- take protective measures to ensure the safety of our receiving and maintenance personnel
- mark the package as appropriate for hazardous materials

APPLICABLE STANDARDS AND DIRECTIVES

We declare herewith that this product complies with the following standards and directives:

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© Heidolph Instruments GmbH & Co. KG