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#### Proper use

The STICKMAX adhesive mat, with dimensions of 20 x 20 cm, makes the use of fixing clips unnecessary in many situations. It is designed for holding glassware (e.g. Erlenmeyer flasks) in place on all types of metal shaking tables and attachments used with shakers and shaker incubators. The glassware used must have a flat, even base. The STICKMAX adhesive mat cannot be used in water bath shakers.

#### Safety instructions

- Always wear personal protective equipment in accordance with the hazard class of the media being worked with.
- The surface of the adhesive mat must be free of grease, dirt, moisture, and oil.
- Do not exceed the maximum load indicated for the device.
- Check all vessels placed on the mat for damage. Never use damaged vessels.
- Ensure that no condensation forms on glassware when used in a shaker incubator.

#### Assembly

The STICKMAX adhesive mat is suitable for use with all types and sizes of shaker tables and shaker attachments (including fixing clip supports). Where necessary, the adhesive mat can be cut to the desired size; guide markings on the underside of the adhesive mat simplify cutting to the right size. Please refer to *Pic 1* for sizing information for IKA shaker tables and shaker attachments.

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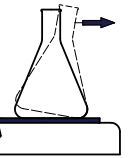
EN

- Degrease the shaker table or shaker attachment (e.g. with isopropanol).
- Moisten the shaker table or shaker attachment with water to allow easy positioning of the adhesive mat.
- Remove the protective film from the side with the writing (underside).
- Submerge the adhesive mat completely in water and then place onto the shaker table or shaker attachment with the "STICKMAX" logo on the underside
- The film of water allows the adhesive mat to be moved around easily for precise positioning. Remove any excess water using a ruler or similar straight edge and mop up with a cloth.
- Allow to dry for 24 hours before first use. Finally, remove the second protective film on the upper side.

#### Use, handling and cleaning

Permissible ambient temperature	C°	5 to 80
Maximum shaker speed	rpm	300
Maximum load	kg	3

Avoid strong or supplemental UV light.

- Clean using a damp, lint-free cloth and a little detergent.
- Disinfect using e.g. 70 percent ethanol.
- To remove the vessel, push or pull gently to the side at the top edge and wait for a few seconds until the vessel slowly releases.
- The adhesive mat can be removed from the shaker table or shaker attachment by pulling to the side diagonally.
- The special properties of the material can be refreshed by placing the mat in water.

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IKA® Works Inc.
IKA® Korea Co LTD